HY-300 Glass bead sterilizers







Product Numbers : HY-300

Product description: HY-300 glass bead sterilizer use a heated glass beads for small lab equipment for fast sterilization products. Sterilizers has compact design and effectively sterilize small solid metal and glassware within ten seconds. When glass beads are heated to about 300 $^{\circ}$ C, it can effectively destroy bacteria and spores. The model is usually used for forceps, scissors, forceps, scalpels, needles, ring vaccination and inoculation needle and commonly used in research labs.

Product Features

- Compact appearance design, digital display temperature at realtime. It can be used in ultra clean workbench.
- 2. Stainless steel inner container, built-in high-performance heating element, high precision of temperature controller and overtemperature protection.
- 3. Inner temperature of container is equipped with high temperature sterilization bead can be heated to about 300 $^{\circ}\text{C}$.
- 4. Small volume, light weight,simple operation, long working life.5 Randomly configurate diameter 1.5 ~ 2.0mm of glass beads.

Technical parameters

Model	HY-300	HY-300D
Temp. Control Range	100°C ~300°C	100°C ~300°C
Temp. Stability	≤±5°C	≤±5°C
Heating Speed	≤25min	≤25min
Heating Well Volume	Ф40 x 80mm	Ф40 x 150mm
Glass Beads Capacity	150g	150g
Cover Material	Stainless Steel	Stainless Steel
Insulation Material	Green Mineral Wool	Green Mineral Wool
Voltage	AC 220 / AC 110V,	AC 220 / AC 110V,
	50/60Hz	50/60Hz
Power	120W	120W
Fuse	250V,2A/3A, Ф 5×20	250V,2A/3A, Φ 5×20
Dimension (WxDxH)	130 x 150 x 200mm	130 x 150 x 240mm
Net Weight	2kgs	2.5kgs

Attention:

- 1, with clean, dry glass beads filled in sterile container. The standard depth below the edge of stainless steel container about 2cm. Please don't use other fillers or liquid contact to container.
- 2. Turn on the power switch and take some time to heat. Then the vessel temperature can reach to the set temperature. Under the conditions of power turning on it will keep the internal temperature.
- 3. Don't touch the glass beads or sterilization container before cooling! Avoid burns!
- 4, Insert the drying equipment needs to be sterilized to sterilized glass beads as much as possible at least 10s.
- 5, Turn off the power switch and the closed the cover after finished to prevent dust and foreign matter.
- 6, Sterilized glass beads must be cleaned at least once per week. Should wait until the instrument cooled completely when getting out the glass beads.
- 7, Reuse after strictly clean and dry. (Get the glass beads out and clean with mild detergent soap, After rinsing ,it can be dried)